

FP4-COL-1550 Installation/Uninstallation Instruction

1. Carefully pick up an FP4-COL-1550 and press it in the opposite side of the fiber connector of any of the fiber adapter models, FP4-FC, FP4-ST, FP4-SC, and FP4-LC. You can use the included screws to hold the part more easily.



2. The rubber gasket on FP4-COL-1550 allows it to stay inside the fiber adapter. Remove the screws if you used them to insert the component in the fiber adapter.



3. Insert the fiber adapter to 818-FA adapter holder. To achieve the most repeatable measurements, use the included screws to fasten the FP4-COL-1550, fiber adapter, and 818-FA.



4. To uninstall FP4-COL-1550, reverse the procedures 1 – 3.